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LED RING ILLUMINATOR WITH POLARIZING FILTER L-120DF

Thank you for purchasing the LED Ring Illuminator With Polarizing Filter L-120DF.
Please read this manual thoroughly before use to insure proper operation and long service.


PACKAGE CONTENTS

※Please check upon receipt.

① Polarizing Filter ... 1

② LED Ring Illuminator ... 1


③ AC Adapter ... 1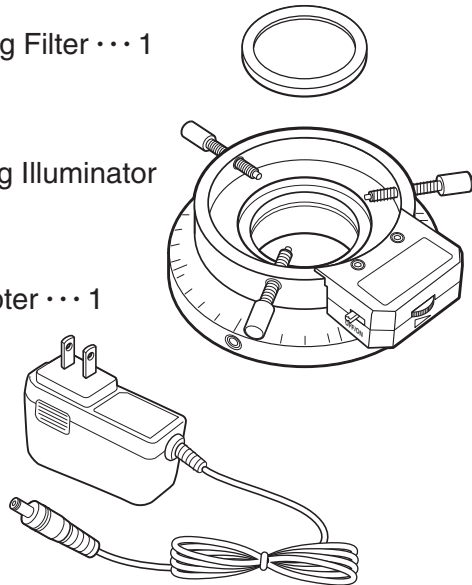


PART IDENTIFICATION

● Polarizing Filter

※Note proper orientation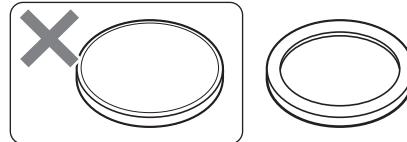

● AC Adapter

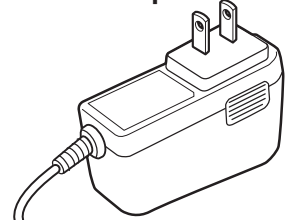● LED Ring Illuminator


Mounting Flange

CLAMP SCREWS
(3 POINTS)

Filter Bezel

※Scale markings are for reference only and do not represent an absolute position or polarization strength.

LED Light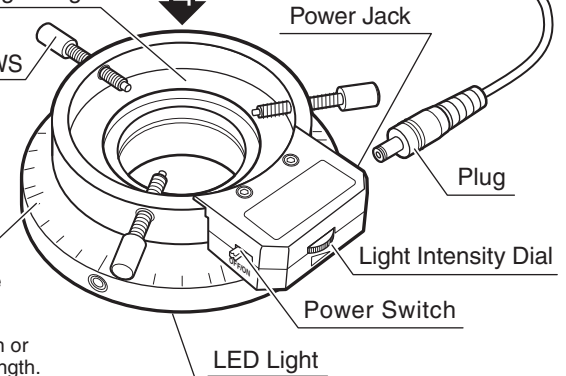

SPECIFICATIONS

Power Supply	AC100-240V 50/60Hz (AC Adapter included)
Power	approx. 5W
Bulb	120x White LED
Color Temp.	approx. 6,500K (daylight)
Illuminance	Max. ~14,000 lx (Measured at distance of 100mm)
Diameter of Microscope Mount	approx. $\phi 30 \sim \phi 60$ mm
Weight	approx. 240g
Dimensions	approx. 90 × 105 × 35mm
Cord Length	approx. 1.4m
Accessories	AC Adapter (Output 24V), Polarizing filter

CAUTION

- Do not use when the unit or LED light is covered with cloth or paper.
- Do not pull or twist the AC Adapter cord. Mishandling of cable may cause failure or result in electrical shock.
- Do not disassemble or modify product; it may cause overheating, fire, or damage to the product.
- Use only as directed; misuse may cause accident or injury.
- Due to product improvements, illustrations shown may vary from actual product.

MOUNTING

※Follow the steps below for proper installation.


- ① Attach the Polarizing Filter to the unit with the proper orientation (Refer to PART IDENTIFICATION above.)
- ② Mount the LED Ring Illuminator to the Microscope Lens Barrel or Adapter provided with the Microscope and secure by uniformly tightening the three Clamp Screws.
- ③ Adjust the polarizing strength by rotating the Polarizing Filter.

■ For models with straight sided Lens Barrel ...

No Adapter is required, just secure directly to Lens Barrel with the Clamp Screws.

■ For models with tapered Lens Barrel ...

Mount the Ring Light Attachment to the bottom of Lens Barrel and secure the Illuminator using the Clamp Screws.

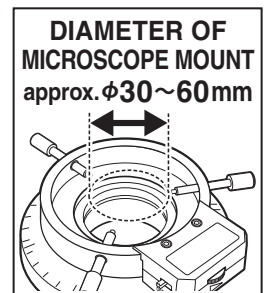Example of mounting:
ZOOM STEREO MICROSCOPE
XTS-SP2